

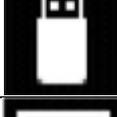
PF7+ and PF7(QC14) ancillary device compatibility guide

Compatibility of attached accessories for PF7+ and PF7 (QC14) are represented using the following icon scheme in the pump HMI:

	Printer	Balance	USB memory stick	Keyboard
Compatibility verified accessories (white icon)				
Compatibility unverified accessories (grey icon)				
Unsupported accessory (red USB icon)				

Verified devices (White Icons)

The following products will be indicated as the corresponding white “verified” compatible accessory icon in the pump HMI (Verified compatible means Flexicon assumes the risk for the accessories compatibility.):

Printer		Epson TM-T88IV Epson TM-T88VI Bixolon SRP-350III
Balance		Mettler Toledo ML-T (104, 204 versions)
USB Memory Stick		Kingston DataTraveler SE9 - 32 GB
Keyboard		Cleanwipe™ Wireless Waterproof Keyboard - SSWKSV099W

Unverified but functioning devices (Grey icons)

Grey icons have an unknown level of compatibility with the PF7+ and PF7 (QC14). All risk of compatibility is assumed by the end user and no promise of compatibility is implied. Any compatibility of unverified accessories is subject to change at any time.

USB flash drive

Product	Comments
SanDisk Ultra Luxe 32GB USB3.1	
SanDisk Ultra Flair 16GB USB3.0	
Samsung Bar plus 32GB USB.31	
Lexar Jumpdrive S75 64GB USB3.1	Must be formatted in FAT32
Corsair Voyager 32 GB USB 3.0	
SanDisk Cruzer Blade 16GB USB 2.0	
Kingston DTIG4/16GB Data Traveler G4 USB 3.1	

Balance

Product	Comments
Mettler Toledo MS-TS	Manual and zero waste batch
Mettler Toledo ME-T	Manual and zero waste batch
Sartorius Practum	Manual filling only

Keyboard

Product	Comments
Logitech MX keys wireless	No issues seen

Non-functioning devices

These may display a red USB device icon and have shown to not function correctly during internal testing.

USB flash drive

Product	Comments
-	Any USB drive that is not formatted in FAT32

Balance

Product	Comments
Mettler Toledo XPR	Dose not work, will not connect to pump
Mettler Toledo XSR	Dose not work, will not connect to pump

Keyboard

Product	Comments
Cleanwipe™ Waterproof Keyboard - SSWKSV099	Wired version of Seal shield keyboard doesn't always initialise properly

Printers

Product	Comments
Epson TM-T88V	Will display a white printer icon but not function correctly